



Agilent 7900 ICP-MS

Specifications and Typical Performance



Raise your expectations with the 7900 ICP-MS

The Agilent 7900 ICP-MS rewrites the rules of ICP-MS. New technologies and a redesigned software platform combine to make the 7900 ICP-MS the world's most powerful, and most automated quadrupole ICP-MS.

Some key components, such as the RF generator and quadrupole mass analyzer, are shared with the field-proven 7700 Series ICP-MS and 8800 ICP-QQQ. However, the Agilent 7900 ICP-MS has been re-engineered from the bench up, with every component optimized for the demands of busy laboratories today, and in the future.



Specifications

| | | |
|---------------------------------------|--|---|
| Sample Introduction System | Peristaltic pump | 10-roller, 3 channels |
| | Nebulizer | MicroMist (borosilicate glass) |
| | Spray chamber | Scott-type double-pass (quartz) Controlled temperature range: -5 °C to +20 °C |
| | Ultra High Matrix Introduction system (UHMI) | Optional |
| Plasma | RF generator | Solid state digital drive 27 MHz Variable-frequency impedance matching 500 W to 1600 W |
| | Torch | One-piece (quartz) 2.5 mm id injector ShieldTorch system |
| | Torch position | Horizontal and vertical position: ±2 mm, in 0.1 mm steps Sampling depth: 3 to 28 mm, in 0.1 mm steps |
| | Mass flow controllers (Ar) | 4: Plasma, Aux., Carrier, Make up/Dilution |
| | 5th gas line for alternative carrier gas | Optional |
| Interface | Sampling cone | 1 mm diameter orifice Standard: Ni-tipped with Cu base Optional: Pt-tipped with Cu base |
| | Skimmer cone | 0.45 mm diameter orifice Standard: Ni Optional: Pt-tipped with Cu base |
| Ion Lens | Lens system | Extraction lens Off-axis Omega lens |
| Octopole Reaction System (ORS) | He (collision) cell gas line | Included |
| | H ₂ (reaction) cell gas line | Optional |
| | 3rd cell gas line (low- or high-flow rate) | Optional |
| Mass Analyzer | Quadrupole | Frequency: 3 MHz Hyperbolic rod profile |
| | Mass range | 2–260 u |
| | Mass resolution | Variable from 0.3 u to 1.0 u |
| | Typical mass calibration stability | < 0.05 u per day < 0.1 u per 6 months |
| | Abundance sensitivity (at Cs) | Low mass side: ≤ 5 × 10 ⁻⁷ High mass side: ≤ 1 × 10 ⁻⁷ |

| | | |
|----------------------|--|---|
| Detector | Configuration | Orthogonal Detector System (ODS) |
| | Detector | Dual-mode discrete dynode electron multiplier |
| | Dynamic range | 11 orders (0.1 cps to 10 Gcps) |
| | Minimum integration time | 100 μ s |
| | Minimum dwell time (TRA mode) | 0.1 ms (no settling time) |
| Vacuum System | Configuration | Three-stage differential vacuum system |
| | Vacuum pump | Single split-flow turbo molecular pump Single external rotary pump |
| | Vacuum pump hose length | 1.5 m, 3 m (optional) |
| Software | Instrument control software | ICP-MS MassHunter Workstation software |
| | User access control software | Optional |
| | Chromatographic software | Optional |
| | Single nanoparticle application module | Optional |
| | Intelligent sequencing software | Optional |
| | Three offline user licenses | Optional |
| | | |

Accessories and Peripherals

| | |
|----------------------------|---|
| Autosamplers | Agilent SPS 4 Autosampler Agilent Integrated Autosampler (I-AS) |
| Sample introduction | Integrated Sample Introduction System 3 PFA Inert Sample Introduction Kit Organic Solvent Introduction Kit Humidifier |
| Speciation kits | LC-ICP-MS Speciation Kits Arsenic Speciation Kit Chromium Speciation Kit Capillary LC Interface Kit GC-ICP-MS Interface |
| Peripherals | Water recirculator Water chiller Optional hood Quiet cover for rotary pump |

Instrument Performance

The factory shipping specifications that are confirmed at the factory represent minimum requirements for shipping approval. The actual performance of the Agilent ICP-MS is invariably much higher. The two tables below provide the typical performance of the Agilent 7900 ICP-MS, together with the factory shipping specifications.

| No Gas mode | | 7900 Factory Specifications ¹ | 7900 Typical Performance ² |
|-------------------------|--------------------------------------|--|---------------------------------------|
| Sensitivity (Mcps/ppm) | ⁷ Li | 55 | 140 |
| | ⁵⁹ Co | | 400 |
| | ⁸⁹ Y | 320 | 600 |
| | ¹¹⁵ In | | 700 |
| | ²⁰⁵ Tl | 250 | 520 |
| | ²³⁸ U | | 720 |
| | Background | <i>m/z</i> 9 | <1 cps |
| Detection limits | ⁹ Be | <0.2 ppt | <0.05 ppt |
| | ¹¹⁵ In | <0.05 ppt | <0.02 ppt |
| | ²⁰⁹ Bi | <0.08 ppt | <0.02 ppt |
| Oxide | CeO/Ce | <1.5% | <1.8% |
| Doubly charged | Ce ²⁺ /Ce | <3% | <2.5% |
| Stability | 20 min | <2.0% RSD | <1.0% RSD |
| | 2 hr | <3.0% RSD | <1.2% RSD |
| Isotope ratio precision | ¹⁰⁷ Ag/ ¹⁰⁹ Ag | <0.1% RSD | <0.1% RSD |

| He Gas mode | | 7900 Typical Performance ² |
|--|-------------------------------------|---------------------------------------|
| Sensitivity (Mcps/ppm) | ⁵⁹ Co | 65 |
| Background | <i>m/z</i> 9 | <0.2 cps |
| Interference reduction factor ³ | ⁵⁹ Co/ ⁵¹ ClO | >30 |
| Oxide | CeO/Ce | <0.5% |
| Detection limits ³ | ⁷⁵ As | <5 ppt |

1. 7900 Factory Shipping Specifications. These specifications are detailed in the Agilent publication: Agilent 7900 ICP-MS, Specifications (Publication number: 5991-3779EN).

2. The typical performance values are not checked during the standard installation.

3. Performed in a matrix of 2% HNO₃ + 0.5% HCl.

Site Requirements and Safety

Dimensions

| | | |
|----------------------------|--------|--|
| Mainframe | Width | 730 mm (main cabinet, excluding peri-pump) |
| | Depth | 600 mm (main cabinet, excluding power cord) |
| | Height | 595 mm (main cabinet, excluding exhaust chimney) |
| | Weight | 100 kg |
| Largest shipping container | Width | 1,020 mm |
| | Depth | 1,120 mm |
| | Height | 1,000 mm |
| | Weight | 148 kg |

Environmental

| | | |
|-----------------------|----------------|-----------------------------|
| Operating temperature | Range | 15–30 °C |
| | Rate of change | <2 °C/hr (max. change 5 °C) |
| Operating humidity | Range | 20-80% (non-condensing) |

Utilities

| | | |
|--------------------|-------------------|---|
| Electricity supply | Voltage | Single Phase, 200-240 V, 50/60 Hz |
| | Current | 30 A |
| Cooling water | Inlet temperature | 15-40 °C |
| | Minimum flow rate | 5 L/min |
| | Inlet pressure | 230-400 kPa |
| Argon gas supply | Minimum purity | 99.99 % |
| | Maximum flow rate | 20 L/min |
| | Supply pressure | 500-700 kPa |
| Cell gas supply | Minimum purity | 99.999% |
| | Maximum flow rate | 12 mL/min for He and 10 mL/min for H ₂ |
| | Supply pressure | 90-130 kPa for He 20-60 kPa for H ₂ |
| Exhaust duct | Vent Type | Single vent, 150 mm diameter |
| | Flow rate | 5-7 m ³ /min |

Regulatory Compliance

| | |
|--------|--|
| Safety | IEC 61010-1:2001 / EN 61010-1:2001 |
| | IEC 61010-2-061:2005 / EN 61010-2-061:2003 |
| | IEC 61010-2-081:2001+A1:2003 / EN 61010-2-081:2002+A1:2003 |
| | Canada: CAN/CSA C22.2 No. 61010-1-04 |
| | Canada: CAN/CSA C22.2 No. 61010-2-061-04 |
| EMC | IEC 61326-1:2012 / EN 61326-1:2013 |
| | Canada: ICES-001:2006 |
| | USA: UL 61010-1 (2nd Edition) |
| ISO | Manufactured at an ISO 9001 and ISO 14001 certified facility |

www.agilent.com

Agilent shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance or use of this material.

Information, descriptions, and specifications in this publication are subject to change without notice.

© Agilent Technologies, Inc. 2015

Published October 22, 2015

Publication number: 5991-6397EN